




## Instructions to Activate an Individual Copy of Mnova.

1. Once you have the license file saved on your disc, run the application on each host. The following dialogs notifying you that a license for the NMR plugin was not found will appear:

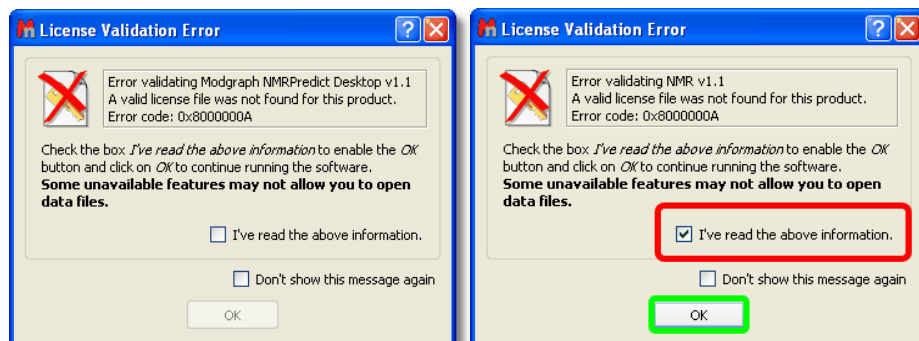

2. Check the 'I've read the above information' boxes, press the 'Ok' buttons and continue running the application. You will obtain the Registration Wizard dialog box, shown below:



3. Click on 'Yes' to display the picture below:





4. Click on 'Activate' (the green square in the picture above), and this will open the 'Specify License File' dialog (see the picture below). Navigate on your Windows browser until you find the license file, and double click on it and select it (see example below). Now you should be able to run the application without receiving an error message dialog.

